

www.precellys.com

NEW GENERATION Tissue homogenizer

SAMPLE PREPARATION & CELL LYSIS



precellys 24
lysis & homogenization

RNA Extraction from Breast Cancer Xenografts and Lymph Node Metastases

at The Institute of Cancer Research

CONTEXT

Within the context of The Institute of Cancer Research, *ex vivo* tumor tissues are being analyzed to study the gene expression of breast tumors and their metastases.

Breast cancer xenografts from different cell lines (MDA-MB-435 and GI-101), as well as their lymph node metastases, were frozen in liquid nitrogen after collection.

MATERIAL

- Precellys®24
- Precellys®24 kit CK14 (small ceramic beads)
- Sample : tumors from breast cancer xenografts and lymph node metastases (frozen)
- Buffer : 600µl of lysis buffer (RTL and β-mercaptoethanol)

PROTOCOL

- Precellys®24 parameters:
5500rpm, 1x20 sec.




RESULTS

In collaboration with The Institute of Cancer Research, McElain Laboratories, Sutton UK.

RNA extraction are performed following 10 different protocols :

Number	Sample	Time	Number	Sample	Time
1	Primary Tumors	10 Sec.	6	Primary Tumors	20 Sec.
2		3 x 10 Sec.	7		Lymph node metastases
3		20 Sec.	8	20 Sec.	
4		20 Sec.	9	20 Sec.	
5		20 Sec.	10	20 Sec.	

RNA gel procedure :
0.5 µg of total RNA diluted in 15µl of denaturing loading buffer (containing urea) denatured and run on a 1% agarose gel in 1x TBE at 140v for 1hour.

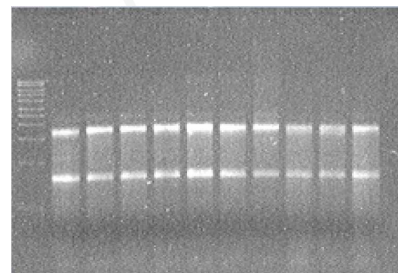

1 2 3 4 5 6 7 8 9 10

Gel electrophoresis analysis after the 10 protocols. Efficiency is validated on the primary tumors and lymph node metastases.

 **The Institute**
of Cancer Research

CONCLUSION

The Precellys® kits allow a quick and effective homogenization of the xenograft tissues, and the total RNA extracted with an appropriate kit following tissue lysis with the Precellys® kits is of good quality.

www.precellys.com
precellys@bertin.fr

Phone : +33(0)139 306 118
Fax : +33(0)139 306 185

bertin
TECHNOLOGIES
mastering innovation

For more details, please contact
precellys@bertin.fr or visit our website

